


2 Step guide to updating drivers for the Digital Microscope on Windows 10.

Step 1: Entering Test Mode

Right click windows start button > Command Prompt (Admin)


In command prompt type this in and enter "bcdedit /set testsigning on"

Restart the computer and it should look like this on the bottom corner.

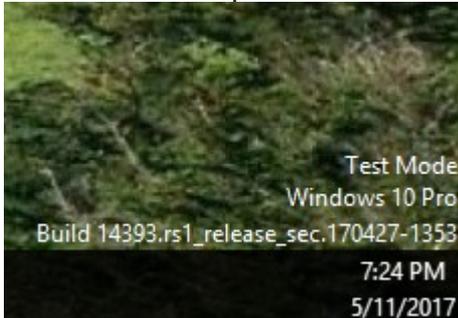



Step 2: Updating drivers

Press Window and search for Device Manager

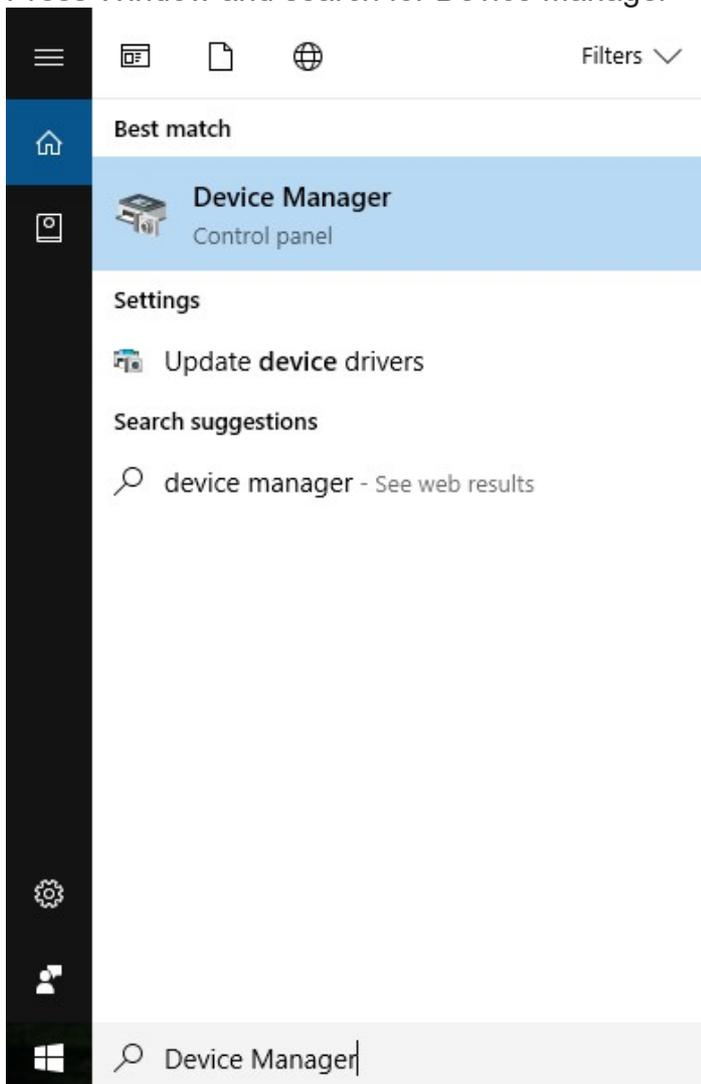

Right click USB2.0 UVC PC Camera and select Update Driver Software

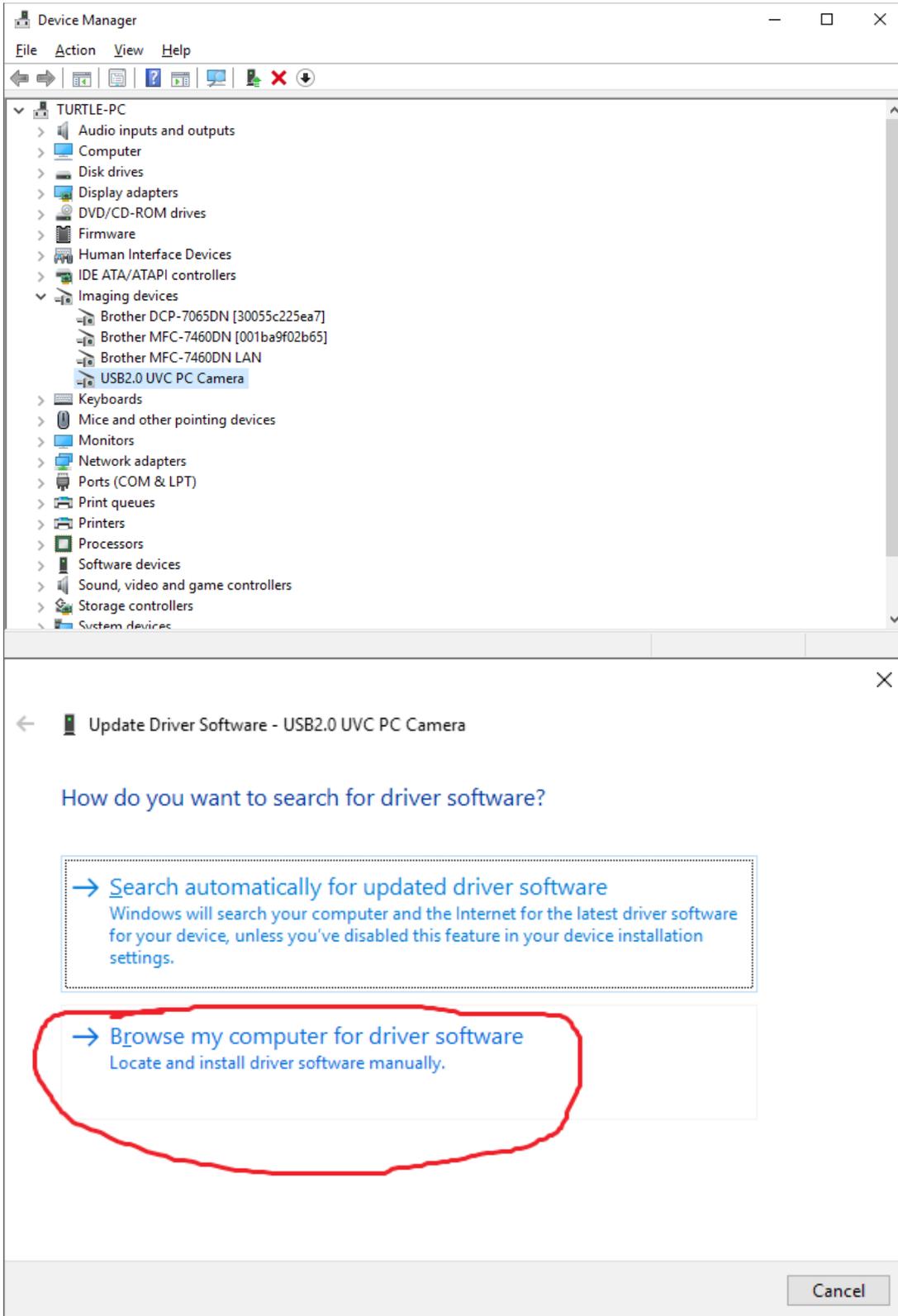

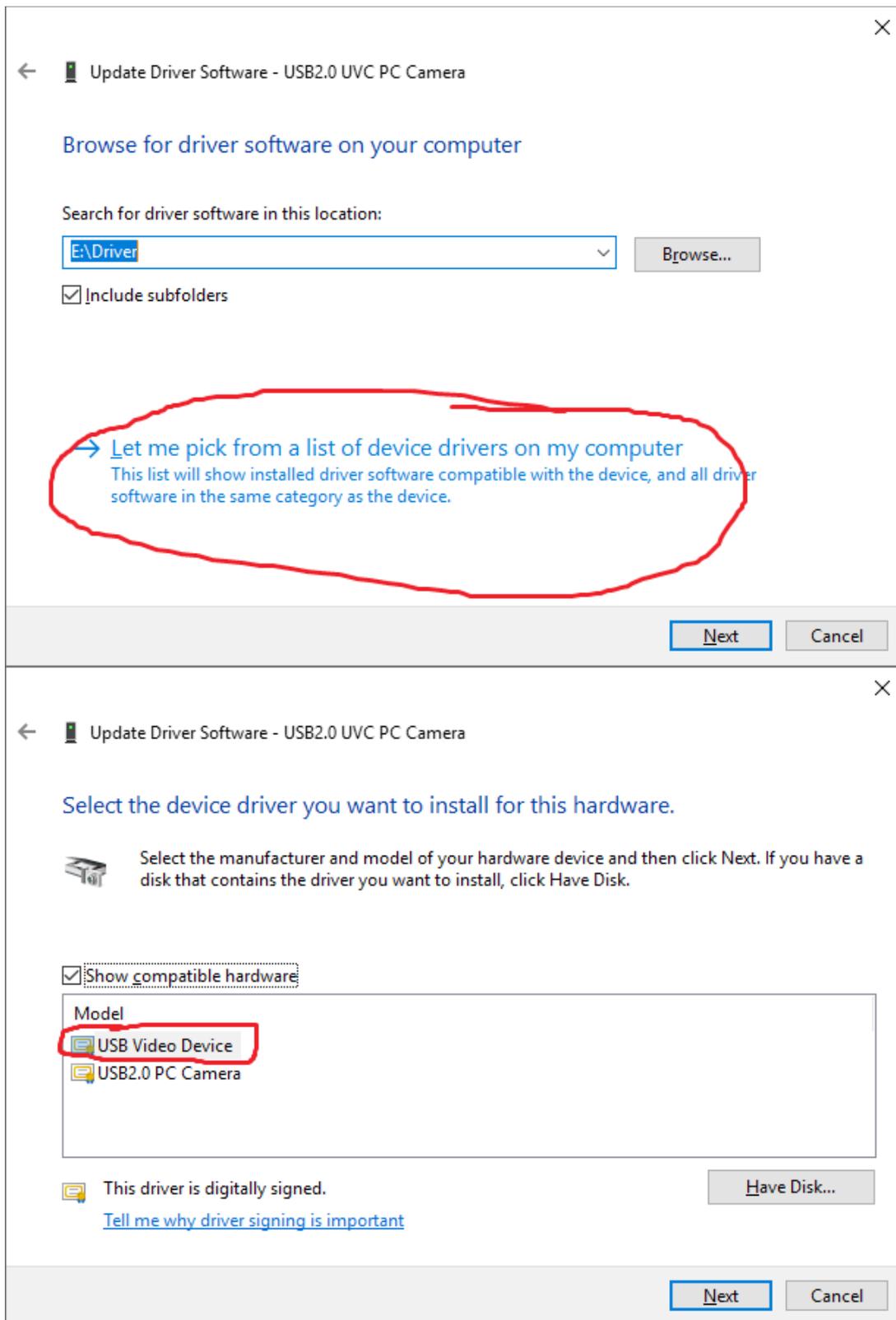

Just 1 mini step left to do!!

Disabling test mode:

Open command prompt again and type in "bcdedit /set testsigning off"

Restart Computer one last time!

Now we are done!